

PI-3 Positive Control

Product No.: 1506738

(For Research Use Only)

■ Product description

SHENTEK® PI-3 Positive Control is intended for use with SHENTEK® Virus DNA & RNA Extraction and Detection Kits, and serves as a positive control to monitor the entire process of PI-3 RNA extraction and RT-PCR detection assay.

■ Kit contents and storage

| Reagent | Quantity | Storage |
|-----------------------|------------------|-------------|
| PI-3 Positive Control | 100 µL × 3 tubes | below -65°C |

PI-3 Positive Control can be stored at appropriate conditions for up to 12 months. Please check the expiration date on the labels.

■ Attentions

1. Thaw one tube completely at 2-8°C or on ice, then vortex and spin briefly.
2. Follow the Positive Control Sample (PCS) preparation protocol outlined in the user guide of SHENTEK® PI-3 Detection Kit and SHENTEK® Bovine Virus Detection Kit (BVDV/REO-3/PI-3/BPV-3/BAV-3).
3. This product is ready for direct use without dilution.
4. Use only after it is equilibrated to room temperature, and avoid repeated freeze-thaw cycles.
5. Do not mix reagents from different batches.
6. Do not use reagents beyond their expiration date.
7. This product is intended for research use only and is not used for clinical diagnosis or treatment. Use with appropriate caution.
8. This product is an armored RNA, which is safe, but it is still recommended to operate it in a biosafety cabinet. Experimental waste generated during its use must be sterilized via autoclaving and disposed of in compliance with medical waste regulations.

■ **Related products**

| Product Name | Product No. | Quantity |
|--|-------------|----------------|
| SHENTEK® Virus DNA & RNA Extraction Kit | 1506730 | 50 Extractions |
| SHENTEK® PI-3 Detection Kit | 1506743 | 50 Reactions |
| SHENTEK® Bovine Virus Detection Kit (BVDV/REO-3/PI-3/BPV-3/BAV-3) | 1506742 | 50 Reactions |

Effective date: 06 Jan. 2025

Support & Contact



Huzhou Shenke Biotechnology Co., Ltd.
www.shentekbio.com
Address: 8th Floor, 6B Building, No.1366 Hongfeng Road, Huzhou 313000, Zhejiang Province, China
E-mail: info@shentekbio.com
Phone: +1 (908) 822-3199 / (+86) 400-878-2189